Declaration of Conformity

EC-Konformitätserklärung
EC-Déclaration de conformité
EC-Dichiarazione di conformità
EC-Declaración de conformidad

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This declaration of conformity is issued under the sole responsibility of the manufacturer.

QUICK GUIDE

TOUCH SCREEN BALANCES

Series: M5

1 Warnings and Safety Instructions

CAUTION!

Please read these instructions for installation and use before starting your work with your new balance. One use for the device other than the one shown in this manual does not guarantee the safety of the product. Keep the instruction manual thoroughly.

The scale complies with guidelines and standards for electrical equipment, electromagnetic compatibility, and the stipulated safety requirements.

If the installation is not performed according to the information provided, or it is made improper use, all warranty rights shall expire.

- **Do not use** the balance in blast risk environment.


- Before starting the unit for the first time, check whether the supply voltage corresponds to the mains voltage.
- To disconnect the unit from the mains power unplug the power cord.


2 Storage

- **Storage temperature:** +5 °C...+40°C
- **Storage humidity** 45% - 75%.
- **Keep package** in the case the balance should be sent back to the factory for service. Disconnect the cables and the accessories for avoiding damages during transportation.
- **Do not place** the balance in extreme temperature and humidity condition and avoid the balance to take strong hit.


2.1 Overview and installation accessories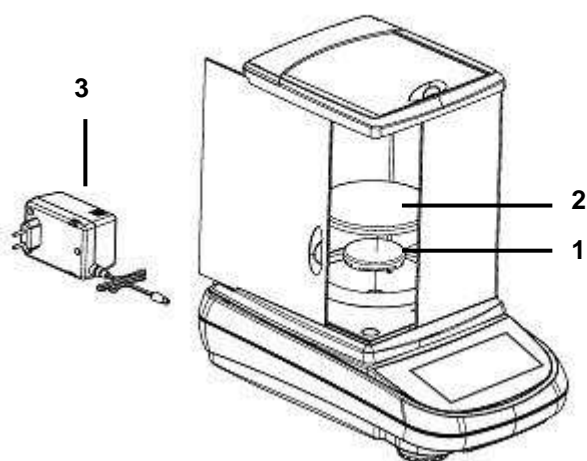


- 1 – Underplate
- 2 – Weighing pan
- 3 – Power supply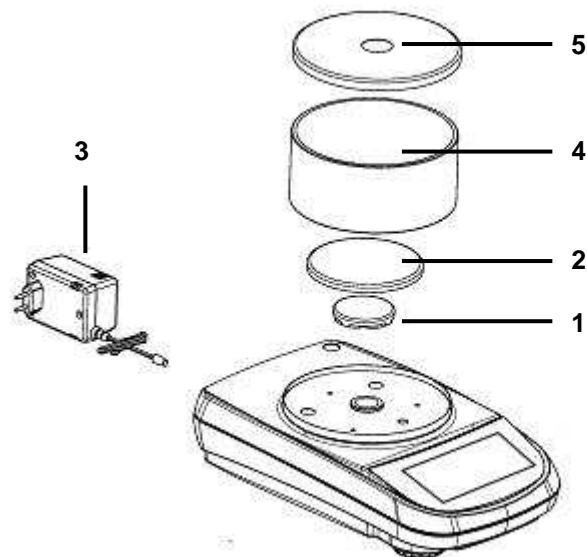

- 1 - Underplate
- 2 – Weighing pan
- 3 – Power supply
- 4 – Cabinet antiventilation
- 5 – Cabinet cover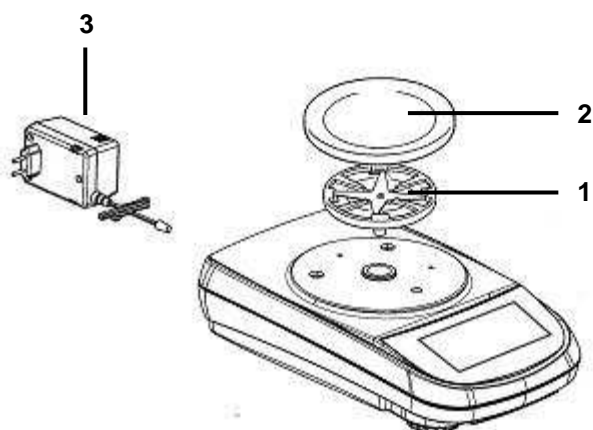



- 1 - Underplate
- 2 – Weighing pan
- 3 – Power supply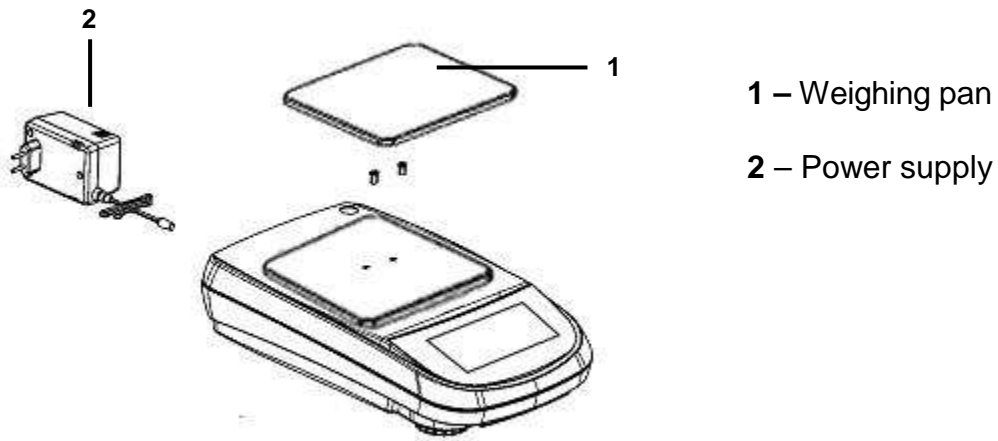

2.2 Place of installation

- **Remove** the balance, calibration weight (if included) and accessories from the cartoon and check for any visible damage of the instrument..
- **Do not install** the balance in a place with air flows, heavy thermal changes and vibrations.
- **Do not use** the balance in blast risk environment
- **The humidity rate** of the balance environment must be between 45% and 75%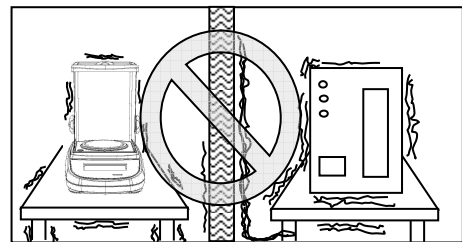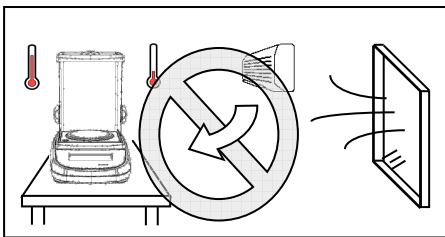

2.3 Initial operation



In models with internal calibration remove the screw in the left lower part of the balance

ATTENZIONE

Level the balance using the level bubble and levelling feet located underneath the case.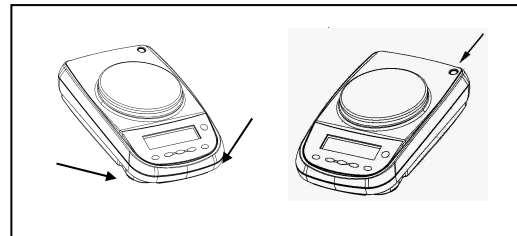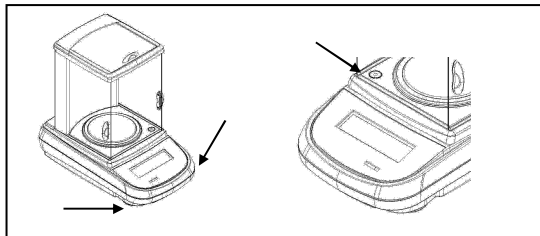

Wait 30 minutes after switching on and calibrate the device after leveling it. For calibration procedures, please refer to the "**Device calibration chapter**". Perform the device calibration whenever it is moved to another place

It is recommended not to drop excess weight objects on the balance weighing plate to prevent damaging it.

Assistance service must be effected by specialized staff and the spare parts used must be original. Therefore, it is necessary to apply to the reseller who sold the equipment.

Plug the jack provided in the connector on the back of the instrument.

Plug power adapter into the power outlet located near the instrument. Do not use non-compliant with current regulations cables / extensions.


N.B. Verify that the power supply indicated on the device nameplate matches the one in use in the country where you are installing the unit.

3 Functionality side keyboard

The balance is equipped with a keyboard placed in the left side part to perform the switching off operations.

3.1 Description

The keyboard has a switching on and off key (1), an operating indicator led (2) and a USB port.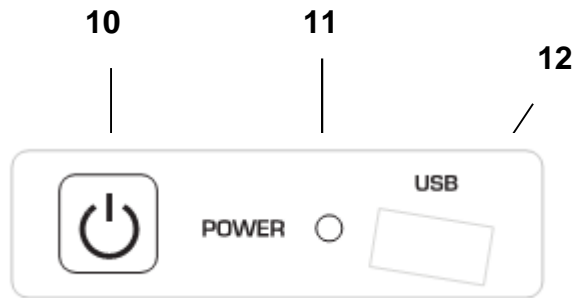

After inserting the power cord, the device will start automatically without having to press the power key.

Use the power key to switch off the device and restart the system when the power cord is permanently attached.

N.B. The phase of switching on and off of the device requires about 20 seconds; during this period, the display will remain switched off.

When powering, wait until you display the home screen before performing other operations.

While in the switching off phase, wait until the green LED turns off before removing the power plug if necessary.


4 Serial interface connections

The device is equipped as standard with a RS 232C interface for connection to a serial printer.

To ensure communication between the device and a printer, you must satisfy the following conditions:


connect the printer via a suitable cable by matching the location of the printer signals with those of the device.

properly set the printer transmission speed and the type of printer according to the printer parameters.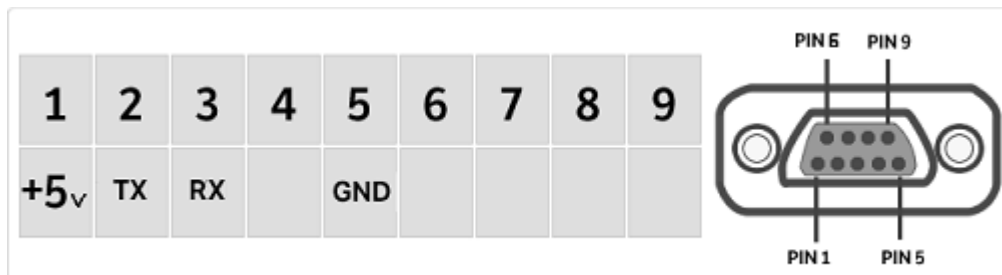

4.1 Display area functions

The balance is equipped with a color display with tap screen functionality that allows tapping on the screen in the active areas to quickly perform the access to the various menus and functionality of the device.

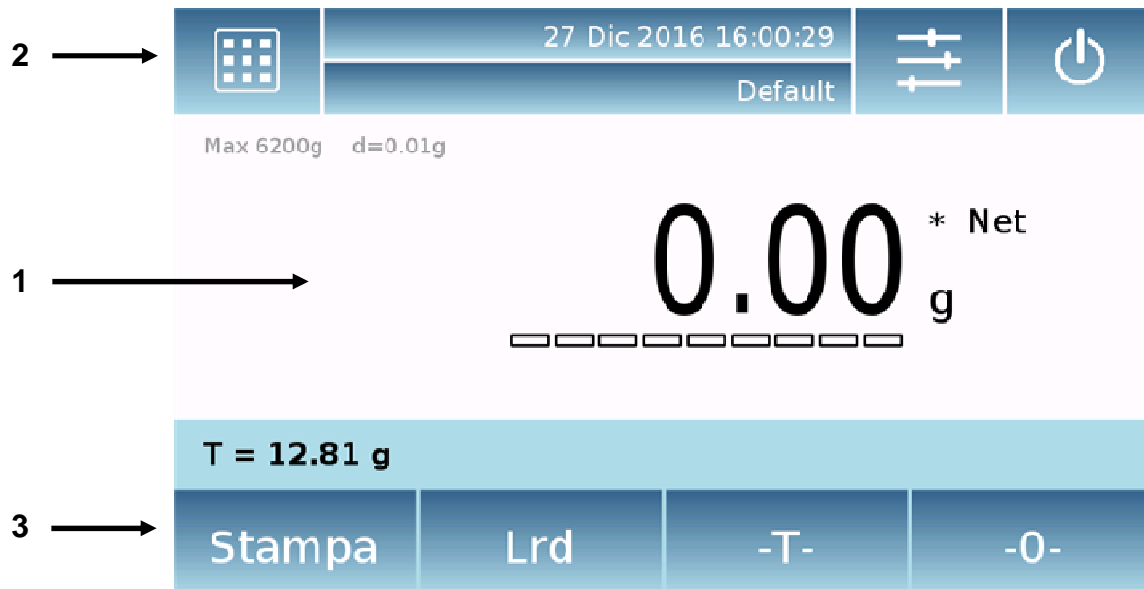1

Area showing the weighing values

g

Display of the measurement unit; tapping the symbol, you can access the menu of the measurement units and select the ones desired.

*

Stability value display.

Net

Indication of the displayed weight. **Net** net value, gross value **Ldr**

T =

Tare value display.

Analog bar showing the magnitude being measured.

Max 6200g d=0.01g

Indication of technical data of the instrument

T = 12.81 g

Information range tare value

2

Bar to display the time and date, to set the weighing parameters, to select the user, to access to drying functions and to switch off the device.

By touching this you access the menu of available functions.

By tapping this area, you access the weighting menu settings, the viewing preferences, the devices and user customization setup.

By tapping this area, you access the switching off menu or enable the standby status of the device,

By tapping this area, you access the user menu to make your selection quickly than desired.

3

Bar to display the function keys.

By tapping the keys displayed in this area, the relative function will be executed. The following are all available keys that may vary depending on the selected applications

Bars of the keys available in simple weighing screens.

Print: key enabling data printing.

-

0-: key enabling the reset key.

-T-: key enabling tare operation.

Lrd/Net: Weight display mode button gross or net.

4.1.1 Instructions for use of the numeric and alphanumeric keypad for data entry

The functionality of the alphanumeric keyboard described below is valid for all of the data insertion operations in the various modes of use of the device.


Mode 1

Mode 2

Mode 3

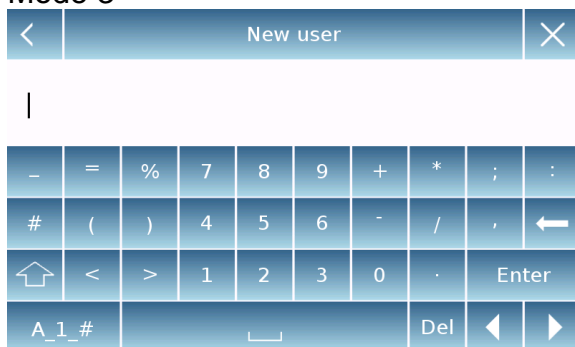Alphanumeric keyboard.

Key to select how to enter the uppercase or lowercase letters

Key to delete the last character entered

Key to delete all the characters entered

Key to move the cursor forward and backward on the string entered

Key to confirm and store the entry

Key to select the keypad characters. By tapping this key, you display randomly the various characters available for the entry.

Mode 1, Mode 2, Mode 3

Key to delete the entry.

<	Tara manuale			×
0 g				
7	8	9	Unità	
4	5	6	C	←
1	2	3	OK	
+/-	0	.		

Numerical keypad

Key to delete the last character entered

Key to confirm and store the entry

Key to delete all the characters entered



Key to delete the entry.

Button for selecting the unit of measurement.

5 Standby, logout, switching off functions

After inserting the power cord into the wall outlet, the device will automatically switch on taking in the basic weighing screen.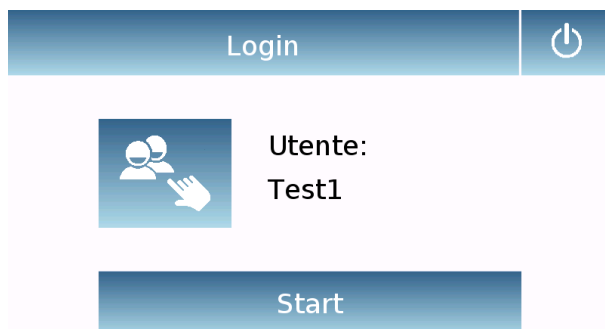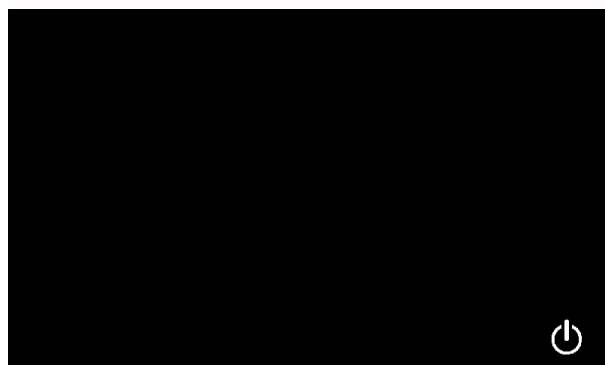

By tapping the area with the symbol  You access the switching off menu.

There are 3 switching off modes:

- Display off.
- Logout.
- Switching off.

Display off

Selecting the display off mode, the device will go in the standby state; the display is turned off, so that only the bottom right command to turn on the system will be enabled.

Tapping the symbol, you will return to the basic weighing screen.

Logout

Selecting the logout mode, you will disable the current user

Tapping "**Start**", it will activate the user and in the case of password protected user it will be asked you to enter it.

Tapping  you can select another user by which you want to switch on the device.

Switching off

Selecting the switching off command, the device will be turned off. To turn it on, you shall press the key on the left of the device or remove the power plug and insert it again.

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